

Mookodi is a low resolution spectrograph and imager, designed to be operated fully robotically. It was designed and built by Liverpool John Moores University (LJMU), and installed on the Lesedi telescope in December 2021.

This (paper) describes the instrument in detail. Please also cite this paper when publishing data collected with this instrument. (for a higher resolution version go [here](#))

Instrument Description

Specifications

- **Camera:** Andor DU934P deep depletion camera, 1024 × 1024 CCD, 13 μm pixels ([CCD performance test sheet](#))
- **Spectral range:** 400.5 – 799.5 nm ^[1]
- **Spectral resolution:**
 - Dispersion:** 3.88 Å/px
 - Resolution:**
 - 80μm (2 arcseconds) slit
 - R = 301.0 @4671.23 Å
 - R = 374.4 @5823.89 Å
 - R = 492.5 @7642.02 Å
 - 160μm (4 arcseconds) slit
 - R = 172 @4671.23 Å
- **Plate scale:** 0.588 arcsec/px
- **Field of view:** 10.07' × 10.07'
- **Slit:** Long slit with narrow (80μm, [2"https://topswiki.sao.ac.za](https://topswiki.sao.ac.za)) and wide (160μm, [4"https://topswiki.sao.ac.za](https://topswiki.sao.ac.za)) sections.
- **Filters:**
 - Filter Slide A:** Full set of SDSS filters (u', g', r', i', z')
 - Filter Slide B:** Longpass order blocking (OB) filter, and clear filter. There are three empty slots in this slide available for future use.

1. [?] Preliminary values based on lab tests at LJMU or original spec. Actual values to be confirmed during commissioning.

Suitable Uses

- Spectroscopy
- Low-cadence imaging: the camera is not a frame transfer camera, and in normal operation exposures are readout after the closure of an iris shutter. The shutter is life-cycle limited and not suitable for high-cadence observations.
- High-cadence imaging: it is possible to deactivate the shutter and take continues short exposure but users should be aware that the camera is not a frame transfer camera and this mode will result in some readout smearing. There is currently also no GPS triggering of the exposures and that the time recorded in the header (Keyword = DATE-OBS, note: is the start of the exposure) is obtained from the CPU time that is NTP synched. However, the header time has been measured to be accurate within at least 100ms of exact time (see [here](#)). **NB. High cadence observing (e.g. multiple continues exposures with exposure times <**

60sec) **SHOULD NOT** be done with the shutter on **AUTO** mode. It is very important to set the shutter to "https://topswiki.sao.ac.za**ALWAYS OPEN**"https://topswiki.sao.ac.za which will deactivate the shutter in the open state but flush the CCD before the start of each new exposure. Failure to do so will likely cause serious damage to the camera's shutter

Data Reduction Pipelines

Bias Correction and Flat-fielding

We now collect automated biases and flats at the start of the night that we collate into monthly master files for users. The files can be obtained [here](#). At that link you can also find a basic reduction python script that can be used to reduce the science files using the appropriate master bias/flat file.

- As an example the script can be used as follow (command arguments should be fairly self explanatory):

```
python mookodi_reduction_using_masters.py --image_dir /home/data/my_target/g --master_bias_file
```

Some important notes:

- In the case of biases, "https://topswiki.sao.ac.za**slow**"https://topswiki.sao.ac.za and "https://topswiki.sao.ac.za**fast**"https://topswiki.sao.ac.za refers to the readout mode and for your science files that can be checked by looking at the "https://topswiki.sao.ac.za**HSHIFT**"https://topswiki.sao.ac.za keyword in the header: slow = 0.05 and fast = 5.0
- Mookodi's derotator cable-wrap does not have a full 360 degree range so the result is that some observations are taken with north up in the images while others flipped 180 degrees with north down in the images. It is roughly correlated to which side of the meridan your observations took place. Your science images can be a mix if a derotation occured during your observation sequence. There are also subtle changes in the telescope throughput for the two orientations so you need to use flats taken at the same orientation as your science images. The supplied reduction script will warn you if you are using flats at a different orientation than your science frames (but still reduce the files). If you are reducing the data yourself with your own software, you can check which orientation your science image was taken at by using the value of the ROTANGLE and ROTSKYPA header keywords. The sky-angle = ROTANGLE-ROTSKYPA. If the sky-angle is close to 0, 360 or -360 your images are with north **DOWN**. If the sky-angle is close to 180 or -180 your images are with north **UP**.

Spectroscopy

The python-based data reduction pipeline **ASPIRED** ([M. C. Lam et. al.](#)) developed for Mookodi's sister-instrument **SPRAT** is compatible and recommended for Mookodi spectroscopic data. General installation and usage instruction can be found [here](#).

More specific Mookodi installation and usage instructions (see included README), with example code and data, can be downloaded [here](#).

Photometry

Differential photometry can be extracted from Mookodi data using a new "https://topswiki.sao.ac.za**SAAO** version"https://topswiki.sao.ac.za, **TEA-Phot-SAAO**, which is a modified **TEA-Phot** pipeline developed by

Dominic Bowman and Daniel Holdsworth ([D. M. Bowman et. al](#))

Installation Instructions

NB. If you use the TEA-Phot-SAAO code to produce results for a scientific publication, we ask that you please adhere to the citation requests of the developers that are on the main repository page (link below).

- These instructions should be identical for a Linux or Mac environment and Python 3.7 or a later version is recommended.
- TEA-Phot-SAAO can be directly downloaded from [here](#) or the repository can quickly be cloned through the following command in the terminal:

```
git clone https://bitbucket.org/saao/tea-phot-saao.git
```

- TEA-Phot-SAAO requires only four additional Python modules. To install (if not already) run the following command in the terminal:

```
pip install numpy matplotlib sep astropy
```

NB. you may need to use "<https://topswiki.sao.ac.za>" if you haven't linked your pip to python 3.6+

Usage Instructions

- To reduce and extract photometry from Mookodi data using TEA-phot-SAAO, in the terminal navigate to the folder that contains the **TEA-Phot.py** file.
- To display basic usage and optional arguments run:

```
python TEA-Phot.py -h
```

NB. you may need to use "<https://topswiki.sao.ac.za>" if you haven't linked your python to python 3.6+

- A typical example to reduce and extract photometry from Mookodi data would be something like this:

```
python TEA-Phot.py SAAO Mookodi MKD --bias MKD --flat MKD --image_dir /full/path/to/science/fits
```

NB. you may need to use "<https://topswiki.sao.ac.za>" if you haven't linked your python to python 3.6+

- Wait for window to appear and select target and comparison star with mouse-clicks.
- Follow further prompts from the pipeline for inputs (aperture size etc.) in the terminal.

Note: TEA-phot will NOT do bias and flat field corrections if not BOTH flat AND bias files are supplied.

Operating Mookodi

Prepare the Lesedi telescope by following the [Lesedi start-up recipes](#). As Mookodi is on the old SHOC port, any relic references to the SHOC port should be taken to mean the Mookodi port.

Mookodi can be operated via a web GUI. Fully scripted observations (incl. auto-focus of the instrument, automatically positioning a target on the slit, and activating auto-guiding) will be available soon.

Operation via the web GUI

Open the web GUI at mookodi.suth.sao.ac.za:5000. The GUI is roughly divided into five sections. From top to bottom: opto-mechanic configuration (lamps, slit, filters etc.), target acquisition (auto acquisition of a user defined sky-coordinate on a user specified x and y pixel position), exposure configuration (exposure time, binning etc.), higher level instrument control (start exposures, auto-focus instrument etc.), and an image display.

1. Opto-Mechanic Configuration:

- ◆ Select the desired configuration for each component. The components moved in/out of the beam via pneumatics (eg. mirrors, slit, grism) and lamps will respond on the push of the button, and indicate when they are busy or in the desired configuration. **To conserve lamp lifetime, do not leave them on unnecessary for extended periods of time.**
- ◆ Select the required filter via the dropdown menu under the Filter A or Filter B headings and press the SET button to activate the filter slides. The box above the dropdown indicates the name of the filter in position for each slide, or BUSY if the slide is still finding its requested position.

2. Target Acquisition: TO DO

3. Exposure Configuration:

Select the desired configuration for each setting using the dropdown menu below the relevant heading, then press the SET button to issue the command to the camera.

- ◆ **Binning:** Options are 1×1, 2×2, 3×3, and 4×4.
- ◆ **Gain:** High (faint objects), medium, and low (bright objects)
- ◆ **Rate (readout rate):** Fast (5 MHz) - read-noise of 13.6 electrons, and slow (0.05 MHz) - read-noise of 3.3 electrons
- ◆ **Image Type:** This has no physical effect on the instrument configuration and only populates a header keyword with the selection (possibly useful for automated data reduction)
- ◆ **Instrument Mode:** This has no physical effect on the instrument configuration and only populates a header keyword with the selection (possibly useful for automated data reduction)
- ◆ **Number of exposures:** The number of exposures to take at the set camera/instrument configuration. **NB. see note on [high cadence imaging in the Suitable uses section at the top of this page](#).**
- ◆ **Exposure time (in ms):** Note the exposure time display does not update when the SET button is pushed, it only updates after the exposure is started. This is a [known issue](#).

4. Higher Level Instrument Control:

- ◆ **Exposure Control:**
 - ◇ **Take exposure/Start:** Starts an exposure with the SET exposure configuration
 - ◇ **Abort:** Aborts the current exposure.
- ◆ **Guider Control:** TODO
- ◆ **Shutter Control:** TODO
- ◆ **Save images:** TODO
- ◆ **Exposure Information:** TODO
- ◆ **Nudge telescope pointing:** TODO

- ◆ **Focus control:** TODO
- ◆ **CCD Temperature Control:** Displays the current CCD temperature, and buttons for cooling down and warming up. The camera takes 10 - 15 minutes to cool, and will stabilise at approximately -70 °C. Cool the camera at the start of observations and warm it up once you are done.

5. Image Display:

- ◆ Contains an image display of the last completed exposure, using [JS9](#).
- ◆ A red and blue circle will appear where the preset values are for the pixel positions for the narrow and wide slit locations. Known bug: only appears after the second image is taken and only displays correctly for 1×1 binning.
- ◆ TIPS: TODO.

Note: It is advisable that DARK frames should be taken with the shutter in the "https://topswiki.sao.ac.zaALWAYS CLOSED" mode. BIAS frames (especially if multiple frames are taken) SHOULD ALWAYS be taken with the shutter in the "https://topswiki.sao.ac.zaALWAYS CLOSED" mode since this will deactivate the iris-shutter in the closed state and prolong the lifetime, see also [high cadence imaging](#) in the Suitable uses section at the top of this page.

Data Access

From a terminal window:

```
scp observer@mookodi.suth.sao.ac.za:/data/lesedi/mkd/yyyy/mmdd/* .<local path to copy to>
```

Password: Same as username shown above with "https://topswiki.sao.ac.zaSao" prefix.

Where yyyy is the year, mm is the month, and dd is the day. The convention is that the data rolls over to the current date at 1200 UTC, hence all the files from a given night are stored in the folder corresponding to the date at the start of the night. e.g. All observations started on the night of 2021/12/11 will be stored in /data/lesedi/mkd/2021/1211/.

Start-up Procedures

It should not ordinarily be necessary for an observer to perform these procedures. However, if a restart is necessary, or if recovering from an unplanned shutdown:

- The Mookodi host machine is set to boot automatically when powered.
- There are three services that need to be run before Mookodi can be operated via the web GUI: opto-mechanics, camera, and web server.

TODO - instructions on how to restart all services in the correct order with a single "https://topswiki.sao.ac.zarestart_services" terminal command

Daily Rounds

This section contains information for the technicians performing daily rounds.

Evening

- Switch on the compressor
- Check that instrument services are running (opto-mech, camera, web service) by opening the web GUI and configuring the instrument as follows:
 - ◆ Slit IN
 - ◆ OB filter SET (Filter Slide B)
 - ◆ 1000 ms exposure SET, Start exposure
- This should produce a noisy image at the camera (visible in the image view window), and all instrument mechanisms should respond as commanded.
- Slit OUT, filter out SET
- If anything fails to respond as expected, follow the troubleshooting procedures.

Morning

- Switch off the compressor
- Check that the camera is warmed up

Troubleshooting

Symptom	Cause	Solution	Who to call
Web GUI does not load	(i) Network connection problem	Check VPN/connection settings	Observer
	(ii) Web server or other services are not running.	Reboot the Mookodi machine and follow the <u>start-up procedures</u> .	IT standby
Filter slide reads 'busy' for excessive amount of time and/or times out to 'N/A'	The filter slide has not been able to register that it has reached its requested position. This has been observed intermittently on Filter Slide A if Mookodi is in an orientation where the filter slide has to extend or retract against gravity.	Command the relevant filter slide to the 'out' position and then back to the required position.	Observer
Pneumatics (cal mirror, slit, grism) do not respond to commands, or read 'N/A' after being commanded	Air supply failure	Check that the instrument air pressure gauge reads between 0.3 - 0.4 kPa. If not, check that the air compressor is on and functioning.	Mechanical/electronics standby

Known Issues

- Exposure time display does not update when the SET button is pushed, it only updates after the exposure is started.
- Filter slide A can time out when trying to reach a commanded position when Mookodi is in certain orientations.