

## Introduction

This page contains additional information for the new SHOC, named "https://topswiki.sao.ac.zaSHOCnWonder" https://topswiki.sao.ac.za, that is now permanently mounted on the 1.9m instrument selector as a secondary instrument. Users should first familiarise themselves with the [Instrument Selector wiki](#).

All the hardware for SHOCnWonder is new, but the underlying control software, GUI, service startup instructions, data folder structures, etc. are identical to the original SHOC systems, so users should also first familiarise themselves with the [SHOC wiki](#).

## Specifications

- **Camera:** [Andor iXon 888 #BV USB3](#) , 1024 × 1024px CCD, 13 μm pixels ([CCD performance test sheet](#))
- **Plate scale:** 0.16 arcsec/px
- **Field of view:** 2.7' × 2.7' (the new system also has a focal reducer so the FoV is roughly double the width of the old system)
- **Filters:**
  - Filter wheel A:** U,B,V,R,I
  - Filter wheel B:** Currently empty (subject to change as these filters are shared with SHOCnAwe and SHOCnDisbelief)

Filter wheel	1	2	3	4	5	6	7	8
A	U	B	V	R	I	--	--	--
B	--	--	--	--	--	--	--	--

## GUI and login details

GUI url (onsite or via VPN access only): <http://shoc74in.suth.sao.ac.za:5000/>

Username and password will be communicated to you.

## Operational changes with SHOCnWonder on the instrument selector

The instrument selector has replaced the old acquisition box, though the guide mirror remains. With SHOCnWonder mounted on the side of the instrument selector, the acquisition procedure is slightly different from when past imaging cameras were mounted on the bottom of the acquisition box. Whereas before, one would centre the guide mirror to allow the beam to pass through the central hole in the guide mirror to the camera, we **no longer move the guide mirror at all**. Instead, a fold mirror is now slid into the beam, covering the hole in the guide mirror that directs the beam to SHOCnWonder. SHOCnWonder is mounted at a 90 degree angle, and is effectively in the "https://topswiki.sao.ac.zainitialise" https://topswiki.sao.ac.za position of the old acquisition & guiding camera. The new guide camera is mounted on linear stages that move in X and Y across the guide star (TCS finderchart) field of view. With the wider field of view, **target acquisition is done using SHOC, not the guide camera**. If the guide camera is centred, it will obscure the SHOC beam, so it should always be positioned away

from centre. The XY slide coordinates at the centre of the field are (7150, 82000).

Quick summary of operational changes:

- Never move the guide mirror
- When using SHOC, keep the fold mirror in the beam at all times
- Do not initialise the XY slides (unless they have encountered a fault and you have first reset them)
- Guide camera is no longer used for acquisition (acquire using SHOC)
- The new guide camera can vignette the science beam - keep it near the edge of the field
- Unlike the older SHOC cameras, there is no option to change the file number manually on SHOCnWonder.

## Observing recipe

### Biases and flats

1. Move the fold mirror into the beam: on the right-hand side of the TCS GUI there is a large green button that will either say "<https://topswiki.sao.ac.za>FOLD MIRROR IS IN BEAM"<https://topswiki.sao.ac.za>, or "<https://topswiki.sao.ac.za>FOLD MIRROR IS OUT OF BEAM"<https://topswiki.sao.ac.za>. If it doesn't say the former, click it to make it so.
2. Park the guide camera in the top-right-hand corner of the guide star field of view (represented by the finderchart on the TCS): click "<https://topswiki.sao.ac.za>GoTo"<https://topswiki.sao.ac.za> in the TCS GUI: XY Slides Control, and enter **X=500, Y=14000**. Wait a few moments for the camera to move then check that it reached the desired coordinates.
3. **For biases:** on the SHOC GUI, **insert the U filter** into the beam. **This is currently necessary due to a light leak** caused by an orange LED, not detected in the U band.
4. Take biases or flats as per the usual SHOC procedures.

### Observing a target

The procedure for acquiring a target with SHOCnWonder on the 1.9m is as follows:

#### On the TCS (tcs74v4), slew to the target:

1. If the guide star finderchart window is open, kill it with the "<https://topswiki.sao.ac.za>"<https://topswiki.sao.ac.za> in the top corner (you need to do this for every target, before a new field can be loaded)
2. Enter your target's RA and Dec into the guide star coordinates GUI, then press ?Set?
3. Press ?TARGET? on TCS GUI
4. Press ?READ? on TCS GUI (RA, Dec will populate from the guide star GUI)
5. Press ?Auto? on TCS GUI (telescope will slew to target)

#### Check fold mirror and focus on TCS GUI:

6. Press so that ?FOLD MIRROR IS IN BEAM?. It should be in beam at all times while using SHOC.

7. If this is the start of your run, set telescope focus (FOCUS F18) to ~2170 (else use the value from last night/previous target)

8. Set Guide "https://topswiki.sao.ac.zaCamera Focus" to ~6000 (enter 6000 in the "https://topswiki.sao.ac.zaGOTO" box and press Enter on the keyboard). Note that the required focus value has been as high as 32000 - further tests are planned to figure out what's happening.

**To ensure that the guide camera is away from the centre of the field -- to avoid obscuring the SHOC beam -- you should select your guide star now:**

9. On guide star finderchart GUI - click on an appropriate guide star in the region outside of the green circle and the red rectangle

•  
Guider field of view with the new instrument selector. SHOCnWonder's field is centred on the green circle. The guide probe enters stage right. Guide stars should be selected from the region outside of both the green circle and the red rectangle.

10. On TCS GUI: XY Slides Control - click on ?Goto GS?

11. On TCS GUI: Exposures & Guiding - select an appropriate exposure time from the "https://topswiki.sao.ac.zaExpose" pull down menu

12. Acquire field on SHOC:

- use SHOC's ?Preview? mode
- adjust telescope pointing
- adjust telescope focus
- adjust guide camera focus if necessary

**Start guiding:**

13. Check that there is a guide star in the Acquisition & Guiding Control image, if not, repeat steps 9 & 10

14. On TCS GUI: Pointer - click ?Guide Star?

15. On TCS GUI: Exposures & Guiding - click ?GUIDE?

**Start photometry:**

16. On SHOC GUI - configure and take science exposures. To cycle through multiple filters, configure SHOC then follow [these instructions](#).

17. To acquire the next target, **do not initialise the XY slides** - this should only be done if there is a fault, in which case the Reset and Initialise buttons will turn red and should be pressed in that order. To move from one target and guide star to the next, simply repeat the procedure above from step 1.

## Data access

There is an automated log being generated as files appear on disk by scraping the fits headers:

[https://www.saao.ac.za/~sbp/74in\\_SHW\\_ObsLogs/](https://www.saao.ac.za/~sbp/74in_SHW_ObsLogs/)

As with the old SHOC systems, data can be copied as follow:

- Open a shell on your local PC (Putty if you're using Windows or a bash/tcsh/sh if on Mac or Linux)
- Copy the data using the following command:

```
rsync -avzP shoc74in@shoc74in.suth.saao.ac.za:"https://topswiki.saao.ac.za/data/74in/shw/
```

- NOTE 1: if the rsync command fails, try it WITHOUT the wildcard (\*) and specify a specific file. Some rsync programs apparently do not support this function.
- NOTE 2: data is also backed-up at the end of the night and on a daily basis to the "https://topswiki.saao.ac.zaastro2015" https://topswiki.saao.ac.za servers so data can also be accessed with the same folder structure from there.

## Scripts for cycling multiple filters

For users that wish to run scripts (e.g. continuously take exposures while cycling a filter sequence), we do not support the old buggy scripting method via the GUI anymore. In fact, on the new system it doesn't work at all so don't even try.

We have developed a new much more stable command line method to run scripts. The instructions are as follow:

- Make sure the SHOC camera is ON and initialised (i.e. you have opened the SHOC web GUI and hit the "https://topswiki.saao.ac.zaTurn Camera On" https://topswiki.saao.ac.za button)
- Setup the camera to your desired readout mode, binning and gain settings (set triggering to "https://topswiki.saao.ac.zaInternal" https://topswiki.saao.ac.za and kinetic series to "https://topswiki.saao.ac.za1" https://topswiki.saao.ac.za)
- ssh to the SHOC PC as the "https://topswiki.saao.ac.zashoc74in" https://topswiki.saao.ac.za user (password is the regular "https://topswiki.saao.ac.zashoc74in" https://topswiki.saao.ac.za user?s password)??

```
ssh shoc74in@shoc74in.suth.saao.ac.za
```

- Navigate to the scripting folder??

```
cd /home/ccd/development/shoc-scripting/
```

- Make a copy of the example script file??

```
cp example_script.txt my_script.txt
```

- Open the newly created script file with nano (or which ever terminal editor you prefer) and edit the script sequence as you desire. The positions you have to enter for the desired filters can be checked on the filter-wheel webgui (click on the blue ?More? text) ??

## SHOCnWonder

```
nano my_script.txt
```

- **NB. very important step** - enter the "https://topswiki.saa.ac.za.shocenv"https://topswiki.saa.ac.za python environment. If you do not do this you will get all sorts of python package import errors when trying to run the script routine??

```
source /home/ccd/.pyenvs/shocenv/bin/activate
```

- Run the script specifying a few fields that will be used to add additional information to the headers (**if your declination is negative, make sure you include the "https://topswiki.saa.ac.za="https://topswiki.saa.ac.za sign and add quotations around it as shown below**):

```
python run_shoc_script.py --observer "https://topswiki.saa.ac.zaJohn Soap"https://topswi
```

- If you get this message on the terminal and you?l also notice the SHOC camera page GUI showing a "https://topswiki.saa.ac.zaError: 500 Internal Server Error"https://topswiki.saa.ac.za:

```
[ERROR] Error running script: TSocket read 0 bytes
```

?\* Then the only way to recover is to restart the SHOC PC and to restart all the services:

```
sudo reboot ?
```

?\* Wait for box the come back online, ssh in again and:

```
restart_services
```

## A comparison between OLD vs. NEW system

Compariing raw unreduced images taken of the same field (07:13:27.974 -15:21:04.21) with similar camera settings for the NEW (left) and OLD (right) SHOC system on the 74-inch.